



Setup Instructions for ER-4500 Energizer/Receivers


If you are adding ER-4500s to an existing installation, shut down VitalView software and unplug all power to the current ER's before adding ER-4500 units.

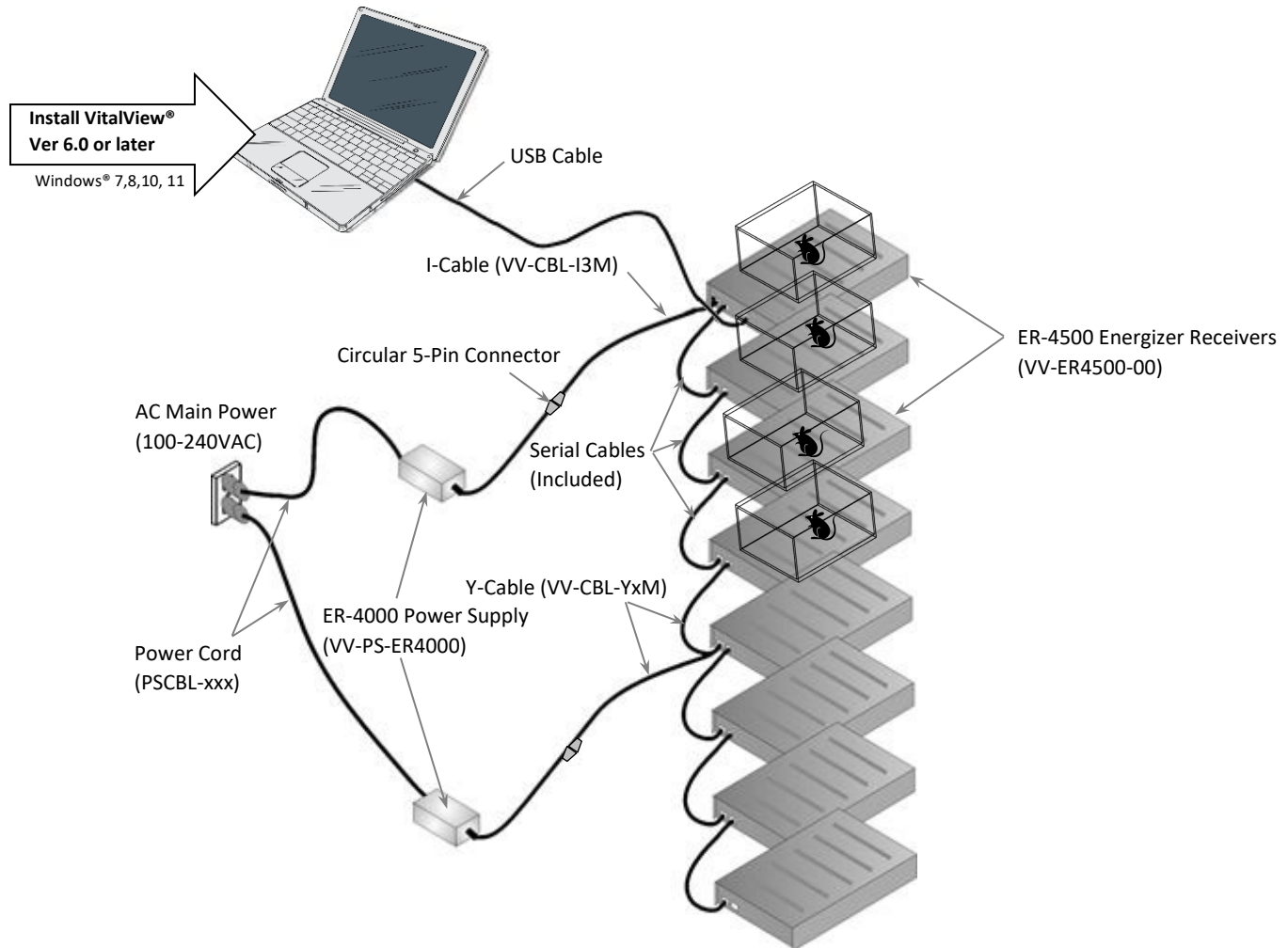

1. Start by placing all ER-4500's and connecting all serial cables.
2. Plug in all power units. When all units have booted up, the display on each will look like the display to the right. Note, all the ID's will be 32.
3. Pick the unit you wish to be 1 and press the square box on the left of the display. This will change the display to look like the one on the right.
4. Press the plus key once and the ID should change to 1.
5. Press the square box on the left to return to the main screen. The ID for this unit is now set to 1.
6. Repeat steps 3 to 5 for each unit using the plus and minus keys to get the ID you wish. **ALL UNITS MUST HAVE A DIFFERENT ID.**
7. You can now install and start VitalView.

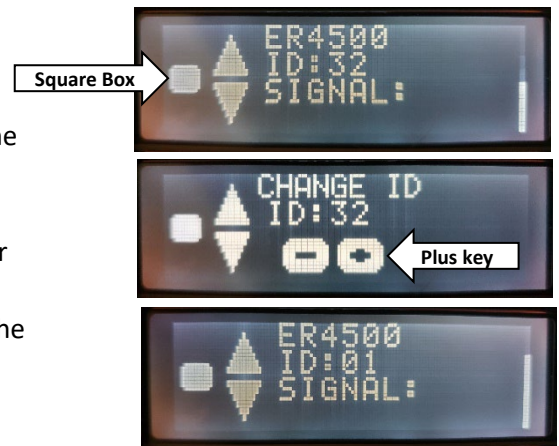

ER4500 Quick Start Guide

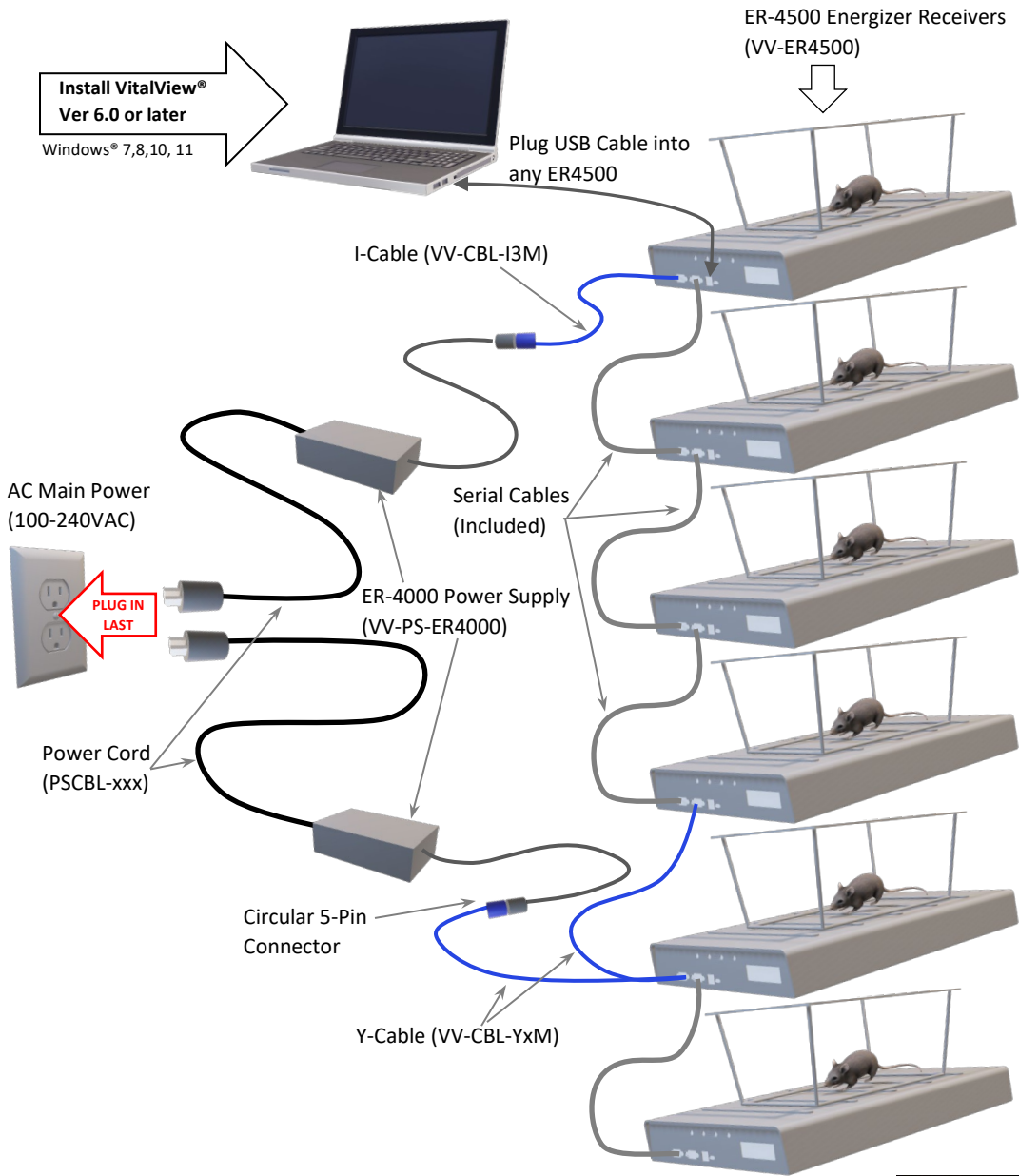

1. Physically arrange ER4500s in desired location in your lab
2. Attach all cabling according to illustration (PLUG IN POWER CORDS LAST)
3. All ER4500s will have default ID 32 (LCD should match Fig 1)
4. You MUST assign each ER4500 a unique ID# before running VitalView
5. E.G. Set ER4500 to ID#1
 - a. Press Enter (■) - LCD should match Fig 2
 - b. Press Plus (+) - ID should change to "01"
 - c. Press Enter (■) - LCD should match Fig 3
6. Repeat step 5 to assign each subsequent ER4500 a unique ID#
7. Run VitalView software when finished



Note: If using ER4500s and ER4000s at the same time, ER4000s must be assigned lower ID#. Minimum spacing between ER4500 and ER4000 is 4 ft (1.2 m). See Vitalview 6 User Manual for detailed setup instructions